

Temperature and level measurement points for the pharmaceutical industry



Negele Messtechnik GmbH

Raiffeisenweg 7 | D-87743 Egg a. d. Günz

Telefon +49 (0) 83 33 / 92 04-0 | Fax +49 (0) 83 33 / 92 04-49

E-Mail: pm@negele.net | www.negele.net

Temperature and level measurement points for the pharmaceutical industry

Negele Messtechnik GmbH has made a name for itself in the food industry as an innovative manufacturer of sensors and process integration systems. In addition, Negele is increasingly establishing itself as an expert in tailored solutions for the pharmaceutical and biotech industries.

In the pharmaceutical industry, a consistently high product quality is critical to the survival of any manufacturer. Apart from reliable and reproducible process monitoring, this requires components that can be cleaned and sterilized with absolutely no residue. The systems and sensors used in production must be designed to reflect this requirement. Negele Messtechnik has brought out two build-in systems for the pharmaceutical industry that it developed specifically to meet these demands.

Aseptic temperature measurement, inline, highly precise and fast – without product contact

Many systems in the pharmaceutical industry exhibit very small nominal diameters. With **PHARMadapt ESP**, Negele offers precisely sized measurement points with nominal diameters from DN 9 for pipe elbows and DN 10 for straight pipe. The exactly defined submersion depth of these sensors ensures measurement precision in both installation situations. PHARMadapt ESP unifies the advantages of a maintenance-friendly, closed system with highly accurate inline temperature measurement and rapid response times. This gives it a significant edge over the commonly used pipe-mounted sensors (Fig. 1).

With a design that is devoid of dead space, the ESP measurement points are easy to clean. In addition, cleaning is supported by the high flow speeds in the CIP process that cause the measurement tip to be actively flushed. The hygienic construction has been awarded certification according to the 3-A 74-03 Sanitary Standard.

The prefabricated build-in pipes can be quickly integrated in the production pipelines in a clean and cost-


effective manner by means of orbital welding. In addition to all nominal diameters up to DN 100 as per the DIN, ISO and ASME pipe standards, the units are also available in different surface qualities (electropolished). All standard certificates in this sector can be provided (including inspection certificate 3.1, surface roughness, and delta ferrite content).

The temperature sensor comes in a variety of models that are compatible with the build-in system. These

Fig. 1: Temperature sensor TFP-58P with build-in system ESP-G DN25



Fig. 2: Temperature sensor TFP-168P with build-in system ESP-G DN10

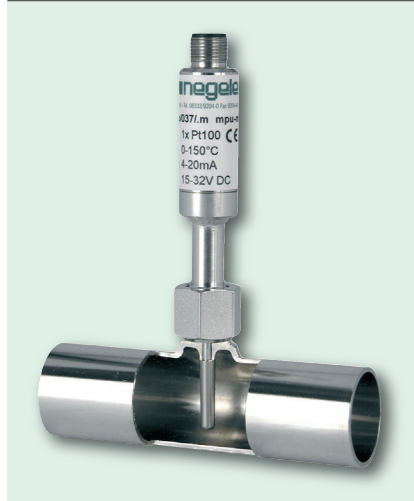


include a standard sensor head for different head transmitters and a miniature version for cramped installation conditions (Fig. 2). A major advantage of the ESP system is the ability to remove the sensor for calibration without opening the process pipelines or detaching the electrical connection. As a result, system maintenance is quick and cost-effective. Another benefit is reliability, since lines cannot be accidentally opened during procedures such as steam sterilization.

Temperature and level measurement points for the pharmaceutical industry

A further build-in system for pharmaceutical applications is the **PHARMadapt EPA** product series with 8 mm and 18 mm holder sizes (Fig. 3). In developing the EPA system, the special requirements and standards of the pharmaceutical industry were taken into account, resulting in a very compact measurement point with an elastomer seal, optimal hygienic design and rapid sensor installation with a clamp ring.

Fig. 3: PHARMadapt EPA build-in system

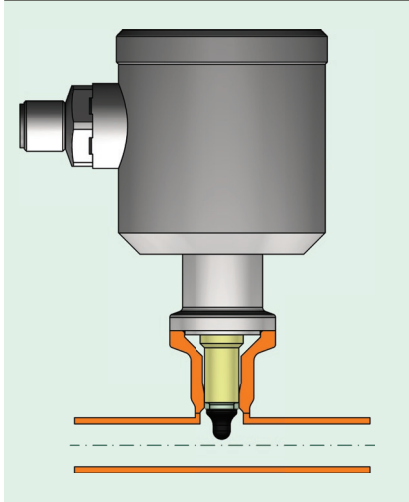

With EPA-8, the user can now not only measure temperatures, but can also record the level in pipes with a nominal diameter of DN 8 and larger. For pipes with diameters larger than DN 25, the EPA-18 installation system is available as a variant with the necessary geometry.

The weld-in pipe and holding sleeves are constructed in such a way that the installation depth of the sensor is always precisely defined. In addition, the configuration is free of dead space and gaps. The measurement point design was developed in

view of the EHEDG guidelines, and the hygienic design of the installation flanges and sensors is certified according to the 3-A 74-03 Sanitary Standard.

Like the ESP build-in system described above, the **PHARMadapt EPA** offers the advantage of prefabricated build-in pipes, enabling the unit to be installed rapidly and cost-effectively by means of orbital welding. Here, too, the DIN, ISO und ASME pipe standards are available with the various surface qualities and certifications.

Temperature measurement with PHARMadapt EPA

The EPA system makes use of temperature sensors that are equipped with weight-optimized and vibration-resistant pharmaceutical sensor heads and that are capable of holding a variety of measurement transducers. These process connections can also hold miniature temperature sensors that feature a M12 plug-in connection or a fixed cable connection (Fig. 4). Components in contact with the product are made of stainless steel 1.4435, for which an inspec-

Fig. 4: Temperature sensor TFP-661 for PHARMadapt EPA



tion certificate 3.1 as per EN 10204 is available. To achieve optimal measuring results, the build-in pipes are available in lengths that match the respective nominal pipe diameters.

Fig. 5: Level switch NCS-6xP for PHARMadapt EPA



Level detection with PHARMadapt EPA

The new capacitive level switches of the NCS series can now be used in these applications (Fig. 5). Designed specifically for the pharmaceutical industry, these devices have been fully reworked and their scope of functions broadened. Thanks to their extended measurement range, the new devices can be used with media that up to now could not be measured. Thus, this new generation of sensors can be used to reliably detect media with a dielectric constant > 2 , such as alcohols and oils. A particularly useful technical innovation is the heated electronics, which prevent the development of condensation in the sensor housing. Other features are a very short response time and the possibility of simulating the sensor states. The new electronics are housed in a weight-optimized sensor head. The standard head design features a window in the cover for viewing the switching state

of the sensor and an M12 plug-in connector with protection class IP 69 K.

The parts of the sensor that are in contact with the process media are made from FDA-compliant materials such as stainless steel 1.4435, PEEK and seal rings of EPDM. The corresponding material verifications are available on request.

With the **PHARMadapt EPA** and **PHARMadapt ESP** build-in systems, Negele Messtechnik GmbH is once again proving itself to be an innova-

tive and flexible provider of measurement point solutions that are a perfect response to customer demands.

Negele Messtechnik GmbH [Hr]

Application example: TFP-58P with ESP-W in a pharmaceutical equipment



Contact



Negele Messtechnik GmbH

Raiffeisenweg 7
D-87743 Egg a. d. Günz

Telephone: +49 (0) 83 33 / 92 04-0
Fax: +49 (0) 83 33 / 92 04-49

E-Mail: pm@negele.net
Internet: www.negele.net

How to find us

